

How to install and set up Repetier-Host (Bioprinting Assisting Software) on a Mac



Overview: The Repetier-Host Software allows Mac users to control their Bioprinters from their computers.













Materials:

You can find the link to download the latest Repetier-Host software [here](#)

Protocol:


1. Navigate to the Repetier-Host page by following the above link or locating the Repetier host link on the <https://cellink.com/support/> download page.

DOWNLOADS

	Category	Version	File
Manual INKREDIBLE	Bioprinter	N/A	Download 
Manual INKREDIBLE+	Bioprinter	N/A	Download 
CELLINK MSDS	Bioink	N/A	Download 
CELLINK A MSDS	Bioink	N/A	Download 
CELLINK Support MSDS	Bioink	N/A	Download 
CELLINK Start MSDS	Bioink	N/A	Download 
Application note	Bioink	N/A	Download 
Python software	Software	N/A	Download 
Slic3r G-code preparation software	Software	N/A	Download 
RAMBo drivers	Software	N/A	Download 
Repetier host	Software	N/A	Download 
INKREDIBLE Software	Software	2.0	Download 

2. Download the version of Repetier-Host for Mac


3. Open the file and drag the icon to your Applications folder as shown in the photo

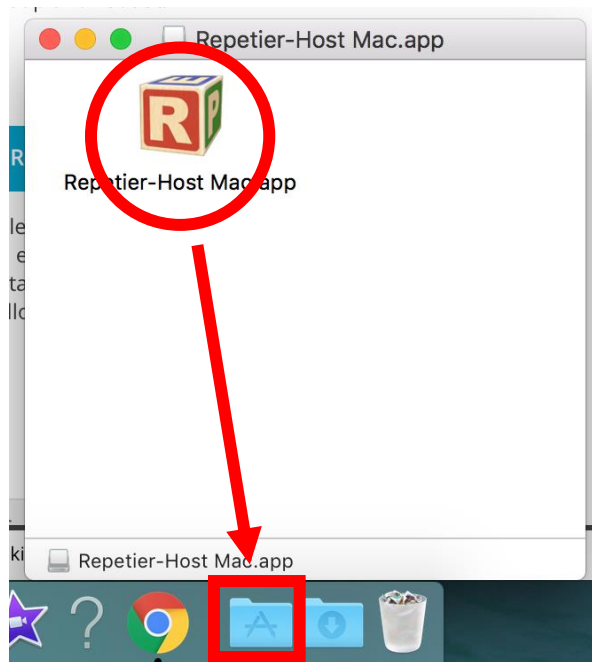


CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00


CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566

4. A security message may appear because the Repetier-Host Application is unfamiliar to the computer. To get around the security settings, open the application from the Applications folder in Finder

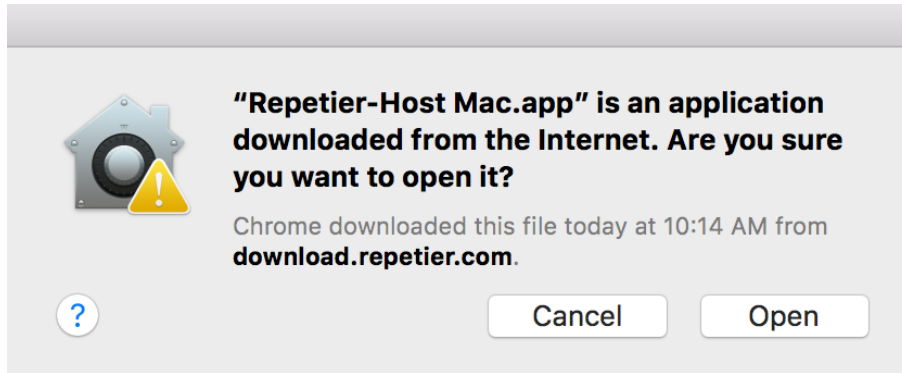


5. Click on "Printer Settings" and ensure that your printer is set to the following parameters

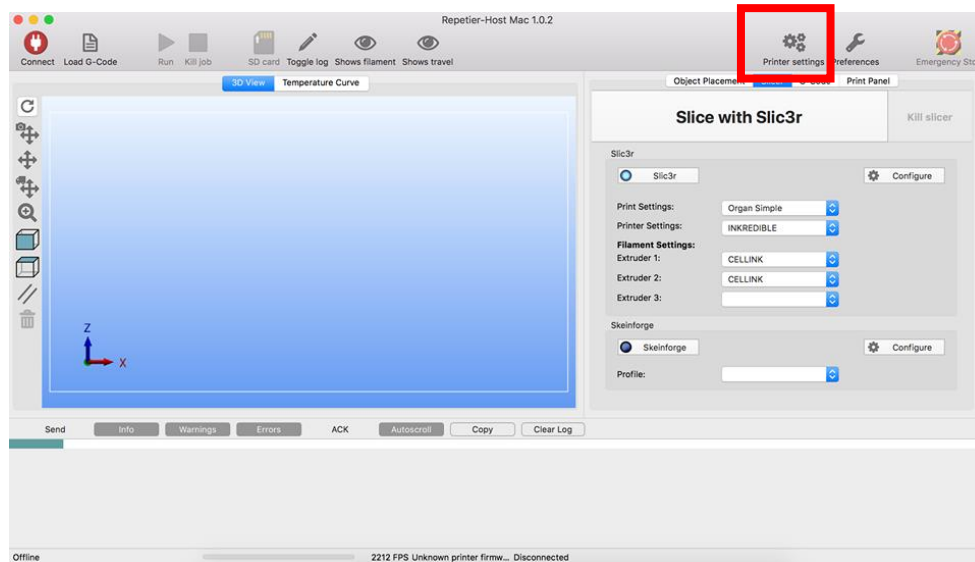

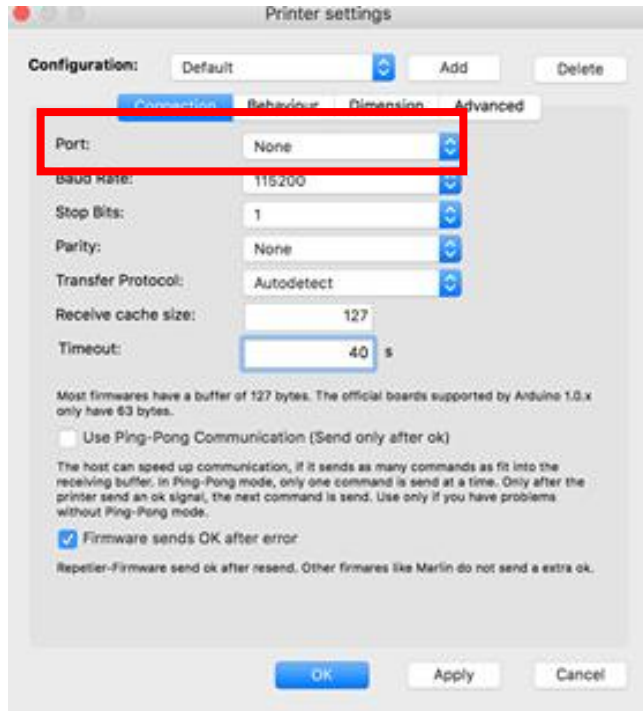


Under port, select the specific port that appears when the computer and printer are hooked up with the USB cable. For example “usbmodem1421”

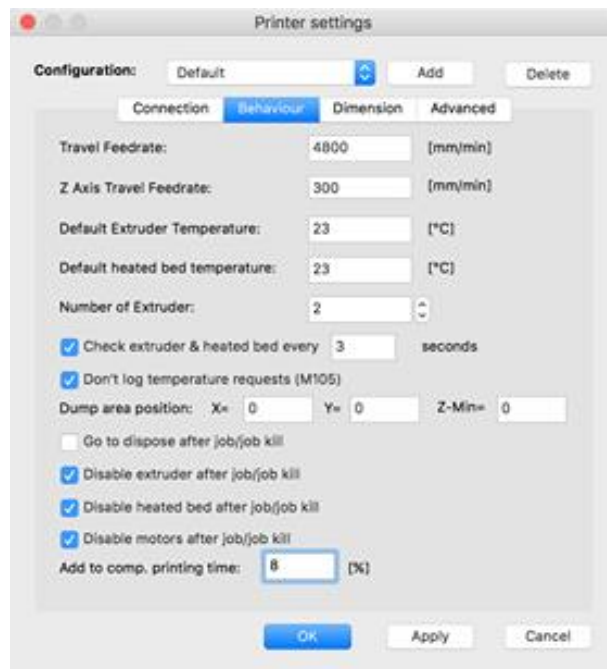

CELLINK AB
Arvid Wallgrens Backe 20
SE 413 46 Gothenburg
Sweden
Phone +46 732 67 00 00

CELLINK LLC
675 W Kendall St.
Cambridge, MA 02142
USA
Phone +1 650 515 5566

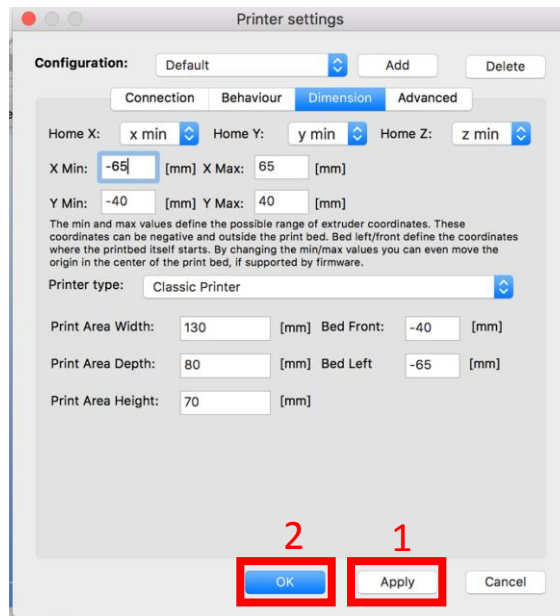

6. Once you are done, press Apply, then click OK

References:

Icon made by Freepik from www.flaticon.com

Icon made by Smashicons from www.flaticon.com